

# IKA

designed for scientists

## IKA R 2850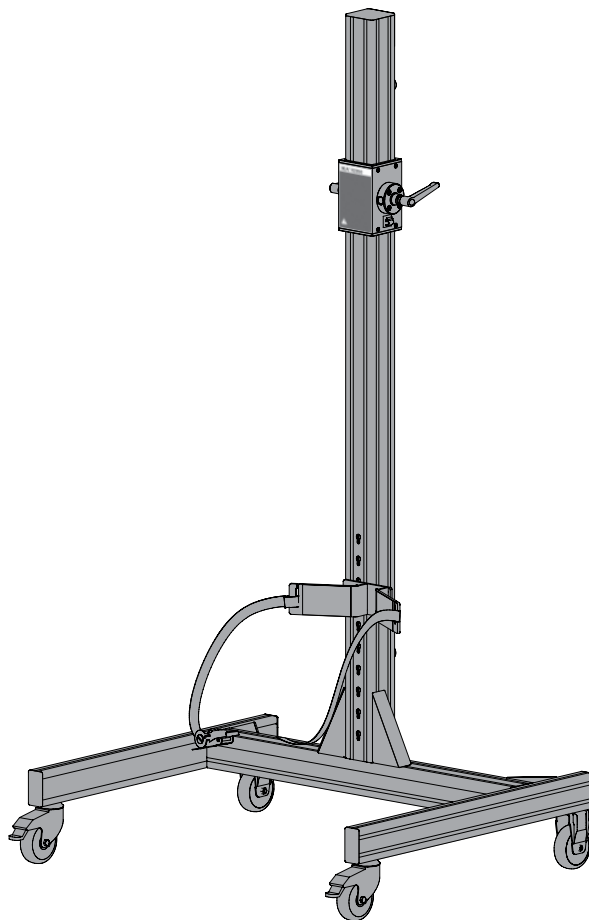

Betriebsanleitung Ursprungssprache	DE	3			
Operating instructions	EN	8	Használati utasítás	HU	75
Mode d'emploi	FR	13	Navodilo za delovanje	SL	80
Руководство по эксплуатации	RU	18	Návod na použitie	SK	85
Instrucciones de manejo	ES	24	Kasutusjuhend	ET	90
Handleiding	NL	30	Lietošanas instrukcija	LV	95
Istruzioni per l'uso	IT	35	Darbo instrukcija	LT	100
Driftsanvisning	SV	40	Ръководство за експлоатация	BG	105
Driftsinstruks	DA	45	Instrucțiuni de utilizare	RO	111
Driftsvejledning	NO	50	Οδηγίες χρήσης	EL	116
Käyttöohje	FI	55	使用说明	ZH	122
Instruções de serviço	PT	60	사용 설명서	KO	127
Instrukcja obsługi	PL	65	取扱説明書	JA	132
Návod k provozu	CS	70			

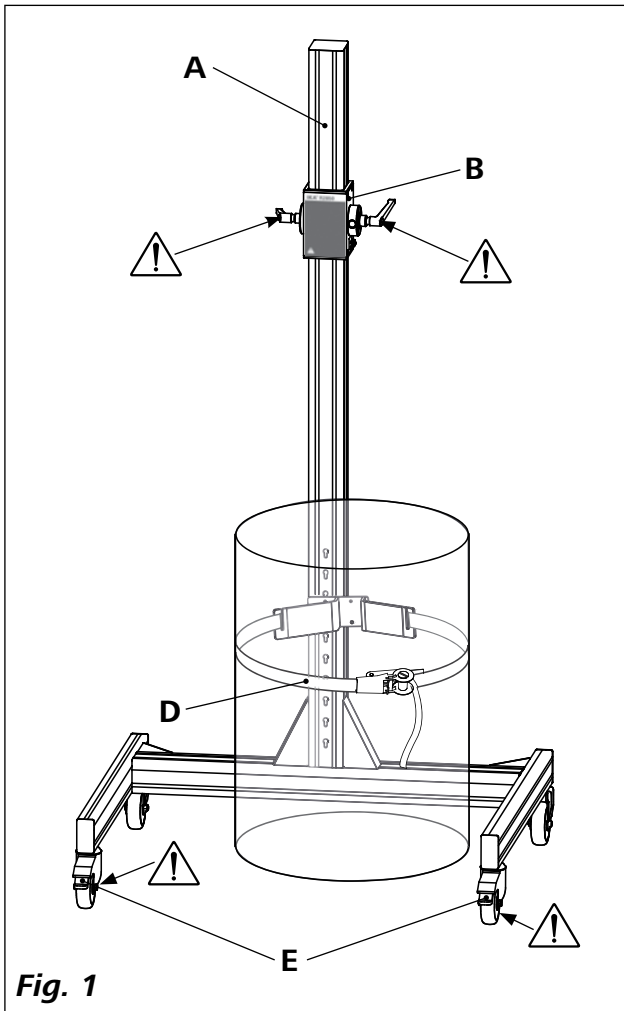**Fig. 1**

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## Contents

	Page
EC-Declaration of conformity	8
Explication of warning symbols	8
Safety instructions	8
Unpacking and transport	9
Intended Use	9
Setting up	10
Alignment	10
Fixing the container	11
Releasing the container lock	11
Maintenance and cleaning	12
Warranty	12
Technical data	12

## EC-Declaration of conformity

We declare under our sole responsibility that the product to which this declaration relates is in conformity with directive 2006/42/EC and the following standard or other normative document: EN ISO 12100.

## Explication of warning symbols

Indicates an imminently hazardous situation, which, if not avoided, will result in death, serious injury.

Indicates a potentially hazardous situation, which, if not avoided, can result in death, serious injury.

Indicates a potentially hazardous situation, which, if not avoided, can result in injury.

Indicates practices which, if not avoided, can result in equipment damage.

## Safety instructions

- **Read the operating instructions completely before starting up and follow the safety instructions.**
- Keep the operating instructions in a place where it can be accessed by everyone.
- Ensure that only trained staff work with the device.
- Follow the safety instructions, guidelines, occupational health, safety and accident prevention regulations.

Please pay attention to the dangerous parts of the equipment in **Fig. 1**.

- Avoid knocks or impact to the stand, stirrer/disperser or accessories.

- Wear your personal protective equipment in accordance with the hazard category of the medium to be processed, otherwise there is a risk of:
- splashing of liquids
  - projectile parts
  - body parts, hair, clothing and jewelry getting caught.

- When used as intended and in accordance with the operating instructions, there is no danger to people, the environment or the device. The manufacturer cannot be held responsible for any damages as a result of incorrect use of the device or non-compliance with the safety instructions.

- Install the stand on a flat, stable, clean, non-slip, dry and fire-resistant surface.
- Prior to each use, always check the stand, stirrer/disperser and accessories for damage. Never use damaged parts.

**WARNING**

Check the mounting position of the stirrer/disperser to avoid the danger of the unit tipping over.

**WARNING**

Observe the maximum permissible load-bearing capacity (see “**Technical data**”).

- You must also follow the safety instructions contained in the operating instructions for the stirrer or disperser.

**WARNING**

Check that the stirrer/disperser is installed correctly before use and at regular intervals:

- the upper and lower end stops should always be tight
- the container should always be held tightly in place using the clamping device
- the lockable castors on the stand must be fixed in place before start-up.

- Because the options for combining products, tools, stirring vessel, trial set-up and medium are nearly endless, user safety cannot be ensured simply with design requirements on the part of the product. For this reason, it may become necessary for users to take other precautionary safety measures. For example, glass apparatus or other stirring vessels that are sensitive to mechanical stress can be damaged or shattered by an imbalance, increasing the speed too quickly or too little distance between the stirring tool and the stirring vessel. Users can sustain serious injury from glass breakage or from the freely-rotating stirrer tool.

**DANGER**

Serious danger of crushing! When adjusting the height of the bracket, ensure:

- your fingers do not get caught in the adjusting slot
- nothing becomes caught between the bracket and devices or other objects attached to the bracket, e.g. the lower end stop or lip of the container.
- When moving the stand, ensure the stirrer/disperser is at the lower end stop position.

**WARNING**

When setting up the stand, ensure that the feet of the unit cannot result in someone tripping over.

## Unpacking and transport

- **Unpacking:**
  - Please unpack the device carefully
  - In the case of any damage a detailed report must be sent immediately (post, rail or forwarder).
- **Transport:**

Once unpacked, the castors can be used to move the stand to its final position.

- **Scope of delivery:**
  - IKA R 2850 Floor stand
  - Tension belt and clip
  - Allen key
  - Operating instructions.

## Intended Use







- **Use:**

The **R 2850** floor stand should only be used to hold, raise and lower stirrers and dispersers. The castors must be locked when the device is in use. The castors must be locked when the device is in use and the container must be held in place using the clips.
- **Field of application:**
  - Laboratories
  - Manufacturing plants



- The safety of the user cannot be ensured:
- if the stand is used in conjunctions with accessories not made or recommended by the manufacturer
  - if the stand is not used for the intended purpose as specified by the manufacturer.

The stand should only be used if it is in perfect working condition. The correct functioning of the stand must be checked at regular intervals.




## Setting up

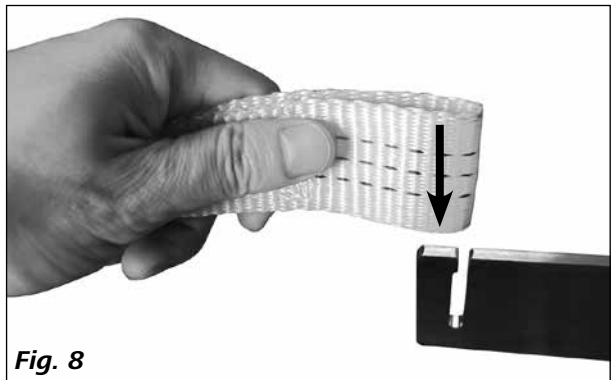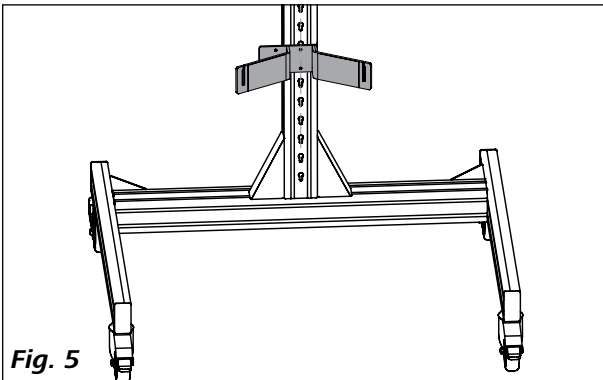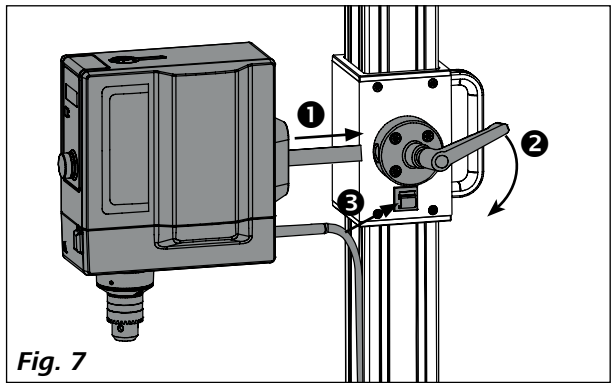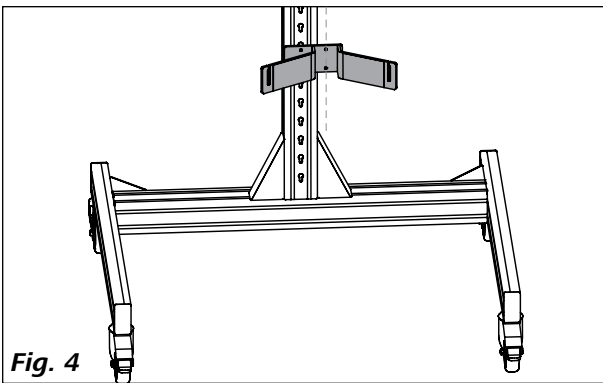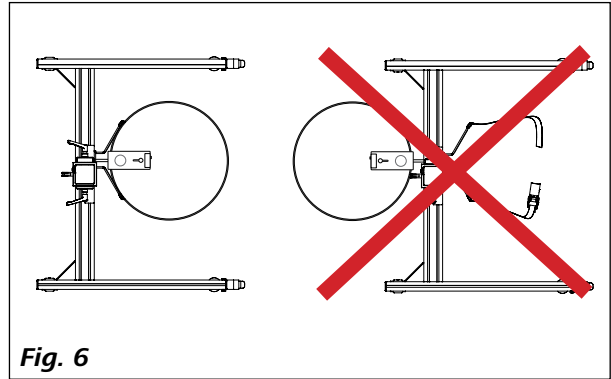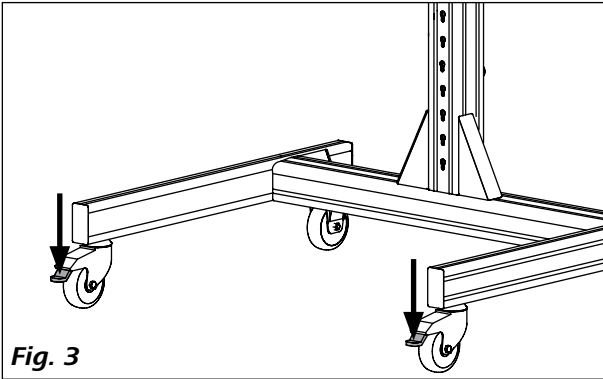

## Alignment
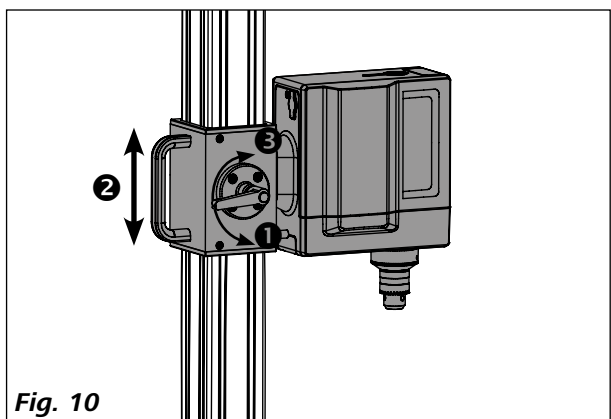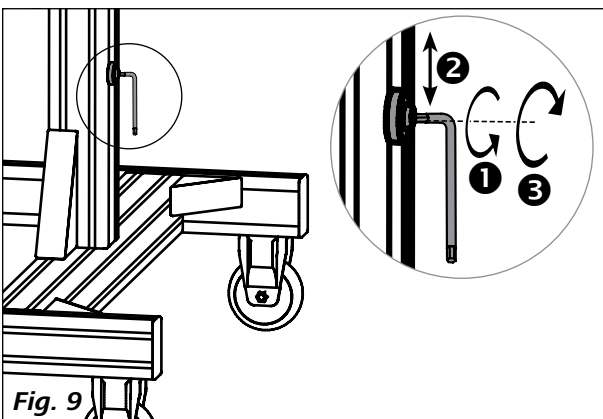
## Fixing the container
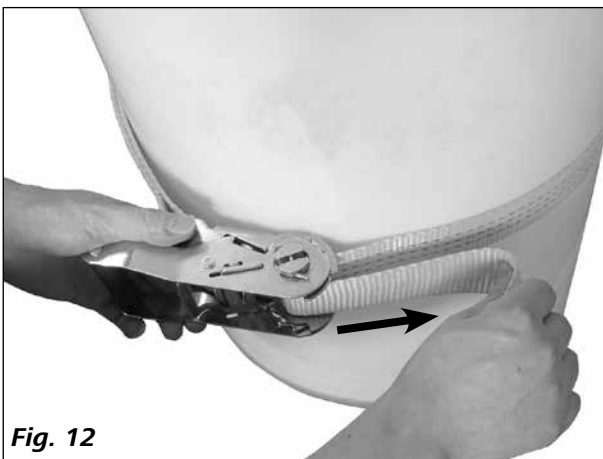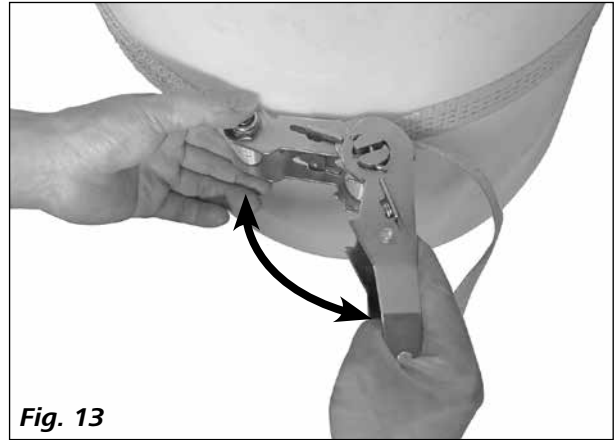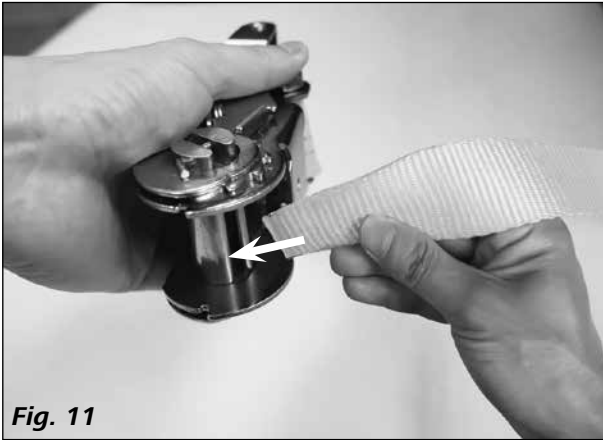
## Releasing the container lock
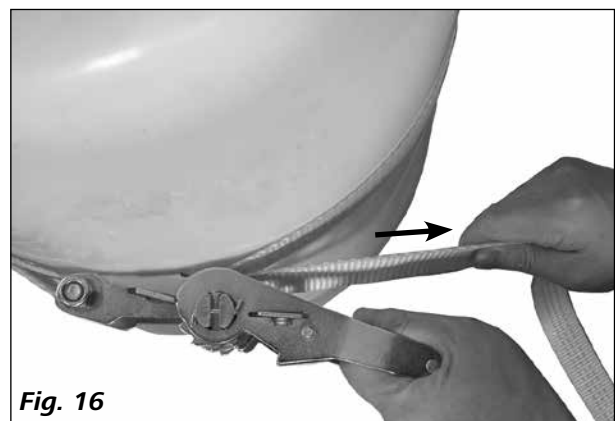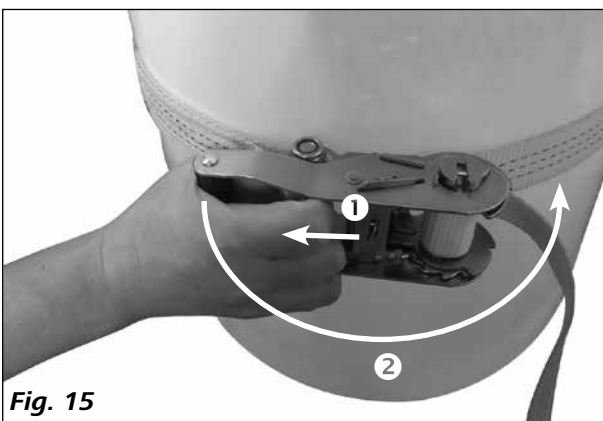
## Maintenance and cleaning

The instrument is maintenance-free. It is only subject to the natural wear and tear of components and their statistical failure rate.

Check regularly that the end stops and brackets are tight.

### **Cleaning:**

Use only cleaning agents which have been approved by **IKA** to clean **IKA** instruments: water containing tenside/ isopropyl alcohol.

Wear protective gloves during cleaning the instruments.

### **Spare parts order:**

When ordering spare parts, please give:

- device type
- manufacturing number, see type plate
- item and designation of the spare part, see **www.ika.com**, spare parts diagram and spare parts list.

### **Repair:**

**Please send in instrument for repair only after it has been cleaned and is free from any materials which may constitute a health hazard.**

For repair, please request the "Decontamination Certificate" form **IKA** or use the download printout of it from **IKA** website: **www.ika.com**.

If you require servicing, return the instrument in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

## Warranty

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

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**IKA India Private Limited**
**IKA England LTD.**
**IKA Vietnam Company Limited**
**IKA Works (Thailand) Co. Ltd.**